

# Nuclear Fast Red Solution

ready-to-use

**42**  
life sciences

WE HAVE THE SOLUTION!

Ready-to-use

Nuclear Fast Red Solution :

- optimized for immunohistochemical (IHC) procedures
- no dilution required prior to use



## Introduction

Nuclear Fast Red Solution is designed for use as a counterstaining solution in immunohistochemistry (IHC) procedures using common chromogens like BCIP/NBT.


Nuclei will be stained red-pink. Cytoplasm will be yellow or unstained.

Specimens counterstained with Nuclear Fast Red Solu-

tion should be dehydrated and mounted with a xylene-based mounting media. This counterstain is not compatible with aqueous mounting media.

## Description

Nuclear Fast Red Solution contains a nuclear fast red stain (a.k.a. Kernechtrot) in stabilizing buffer. No dilution is required prior to use.

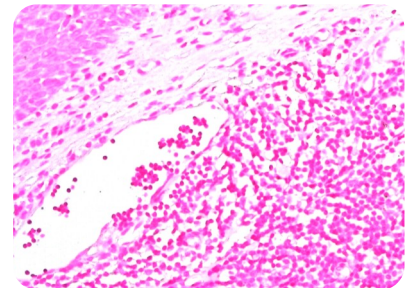

Counterstain with *Nuclear Fast Red Solution* on tonsil section.

## Order Information

Prod. No.	Product
S-99031-1000	Nuclear Fast Red Solution - 1000 ml

42 life sciences GmbH & Co. KG

Fischkai 1  
D-27572 Bremerhaven  
Germany  
www.42ls.com  
info@42ls.com